


Instructions for installing WOLFViewer in Personal Computers

WOLFViewer software can be installed in personal computers for data analysis. However, WOLF instrument operation can only be performed with the computer provided by NanoCollect with the system.

1. To install the WOLFViewer software in a personal computer, click on the following [link](#).
2. Within the link, there will be a .zip folder downloaded. This .zip folder will contain a 'WOLFViewer' folder and 1 Java.exe file. This .exe file needs to be installed on your Windows system for the WOLFViewer software to run properly.
3. Once the files are downloaded, execute the jdk 8U221 Windows X64.exe file.
4. Copy+Paste or drag the WOLFViewer folder (containing the WOLFViewer.exe file and all necessary log files) from the downloads folder to the desktop of the destination computer.
5. Lastly, launch the WOLFViewer software by opening the WOLFViewer folder and double-clicking on the WOLFViewer application.

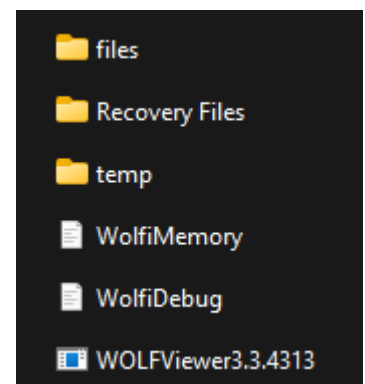

NOTE: The WOLFViewer software will take about 2 to 3 minutes to launch (Figure 2). A message stating "software is not connected, please click connect in the Connection Status panel to reconnect" will be prompted. This is normal as the PC will not be connected to a WOLF instrument. Select Ok. The WOLFViewer launching window will close and the software will open about 10 seconds later.

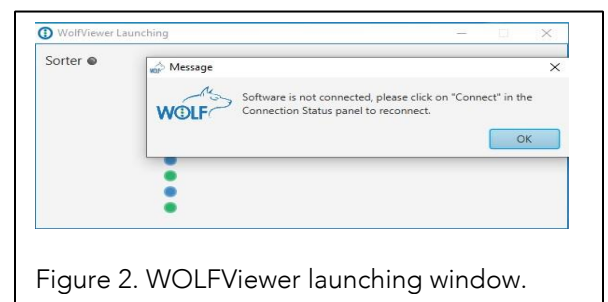


Figure 2. WOLFViewer launching window.

Once the software opens, load your FCS files and enjoy the analysis!

If you need any assistance during this process, please contact Technical Support at (877) 745-7678 Ext. 3 (Office Hours: 6:00AM – 5:00PM Pacific Time).